

Refractometers

Digital Refractometer KERN ORM



Transport and storage case



Rear view, screw-on battery compartment cover

Digital measurement of refraction index for universal application

Features

- The KERN ORM refractometers are accurate and universal maintenance free digital handheld refractometers
- They are characterized by their easy-using and robustness
- The typical and practical design is suitable for a quick and convenient everyday use
- The large, easy-to-read display with integrated temperature display supports the user to reliably determine the measurement
- The integrated automatic temperature compensation (ATC), avoids the manual conversion of the measurement. This allows a quick and efficient usage of the instrument
- Rapid, user-friendly calibration of the refractometer is possible at any time using standard commercial distilled water
- The refractometers from the KERN ORM range are protected to international IP65 protection class, against dust and water splashes. After use, you can rinse the refractometer under running water
- Mean value measurements possible
- The following accessory-parts are included:
 - Prism cover lid
 - Pipette
 - Storage box
 - 1 × AAA battery
 - Screwdriver

Technical data


- Measurement temperature: 0 °C – 40 °C
- Overall dimensions W×D×H 121×58×25 mm
- Net weight approx. 289 g
- Power supply: 1 × AAA (1,5 V)
- Lifetime of the battery:
approx. 10.000 measurements
- ATC (Automatic Temperature Compensation)
- Minimum sample volume: 4 drops
- Automatic energy management
(AUTO-OFF after 60 seconds)
- Mean value measurement (15 measurements)

Accessories

- Calibration liquid




Also available with calibration certificate, see page 108!

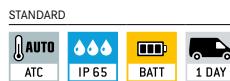

Digital Refractometer KERN ORM-B · ORM-R · ORM-SU

Scope of application: Basic measurements for Brix and refractive index

The following models are particularly suitable for basic measurement where the result is required in Brix or refractive index. They are used to determine the sugar content in food or for monitoring processes in the industry (coolant monitoring, water-based mixtures). Alternatively the display can be switched to show Brix or the refractive index.

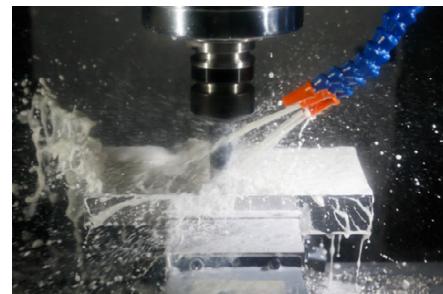

The main scope of applications is:

- Industry: Monitoring of lubricants in machines and quality control
- Food industry: Beverages, fruits and sweets
- Agriculture: Determination of the degree of ripeness of fruit for quality control in harvesting
- Restaurants and large-scale catering establishment



Model	Scales	Measuring range	Accuracy	Division
KERN				
ORM 50BM	Brix Refractive index	0 – 50 % 1,3330 – 1,4200 nD	± 0,2 % ± 0,0003 nD	0,1 % 0,0001 nD
ORM 1RS	Brix Refractive index	0 – 90 % 1,3330 – 1,5177 nD	± 0,2 % ± 0,0003 nD	0,1 % 0,0001 nD

Scope of application: Sugar

The following models are particularly suitable for direct measurement of different types of sugar. These are used to determine the content of the respective type of sugar in water-based liquids. It is possible to switch between the four different scales.



The main scope of applications is:

- Food industry: Beverages, fruits and sweets
- Agriculture: Determination of the degree of ripeness of fruit for quality control in harvesting
- Restaurants and large-scale catering establishment



Model	Scales	Measuring range	Accuracy	Division
KERN				
ORM 1SU	Fructose	0 – 69 %	± 0,2 %	0,1 %
	Glucose	0 – 60 %	± 0,2 %	0,1 %
	Brix	0 – 90 %	± 0,2 %	0,1 %
	Refractive index	1,3330 – 1,5177 nD	± 0,0003 nD	0,0001 nD
ORM 2SU	Lactose	0 – 17 %	± 0,2 %	0,1 %
	Maltose	0 – 16 %	± 0,2 %	0,1 %
	Dextran	0 – 11 %	± 0,2 %	0,1 %
	Brix	0 – 50 %	± 0,2 %	0,1 %

Digital Refractometer KERN ORM-HO · ORM-NA · ORM-SW

Scope of application: Honey

The following model is particularly suitable for the measurement of the water content in honey according to the International Honey Commission (IHC2002) and "degrees Baumé" to determine the relative density of liquids. Alternatively the display can be switched to show Brix or the refractive index.



The main scope of applications is:

- Beekeeping
- Honey production

Model	Scales	Measuring range	Accuracy	Division
KERN				
ORM 1HO	Brix	5 – 38 %	± 0,2 %	0,1 %
	Baumé	33 – 48 °Bé	± 0,2 °Bé	0,1 °Bé
	Water content	0 – 90 %	± 0,2 %	0,1 %
	Refractive index	1,3330 – 1,5177 nD	± 0,0003 nD	0,0001 nD



Scope of application: Salt

The following models are particularly suitable to determin the concentration of NaCl (salt) in water and seawater. This is often used for the preparation and for the cooking of sauces, bases for pastries, the production of brines (e.g. for white cheese) and the preparation of seafood and marinades for meat. Alternatively the display can be switched to show Brix or the refractive index.



The main scope of applications is:

- Food industry
- Restaurants, and large-scale catering establishment, canteens
- Fisch farm

Model	Scales	Measuring range	Accuracy	Division
KERN				
ORM 1NA	Salt content (NaCl) %	0 – 28 %	± 0,2 %	0,1 %
	Salt content (NaCl) ‰	0 – 280 ‰	± 2 ‰	1 ‰
	Salt content (specific gravity)	1,000 – 1,220	± 0,002	0,001
	Brix	0 – 50 %	± 0,2 %	0,1 %
	Refractive index	1,3330 – 1,4200 nD	± 0,0003 nD	0,0001 nD
ORM 1SW	Salt content seawater	0 – 100 ‰	± 2 ‰	1 ‰
	Chlorine content seawater	0 – 57 ‰	± 2 ‰	1 ‰
	Salt content (specific gravity)	1,000 – 1,070	± 0,002	0,001
	Brix	0 – 50 %	± 0,2 %	0,1 %
	Refractive index	1,3330 – 1,4200 nD	± 0,0003 nD	0,0001 nD



Digital Refractometer KERN ORM-AL · ORM-BR · ORM-WN

Scope of application: Beer/alcohol

The following models are particularly suitable for determining the sugar content of the original wort of beer in its unfermented state. The value can be read straightaway, without having to be converted, using the Original gravity (specific weight) and Degrees Plato scales. In addition, the percent by volume and percent by mass scales can be used to determine the alcohol content of clear spirits.

The main scope of applications is:

- Beer brewers
- Alcohol production



Model	Scales	Measuring range	Accuracy	Division
KERN				
ORM 1AL	Percentage by mass	0 - 72 %	± 1 %	1 %
	Percentage by volume	0 - 80 %	± 1 %	1 %
	Brix	0 - 50 %	± 0,2 %	0,1 %
	Refractive index	1,3330 - 1,4200 nD	± 0,0003 nD	0,0001 nD
ORM 1BR	Plato	0 - 31 °P	± 0,3 °P	0,1 °P
	Original gravity (specific weight)	1,000 - 1,130	± 0,002	0,001
	Brix	0 - 50 %	± 0,2 %	0,1 %
	Refractive index	1,3330 - 1,4200 nD	± 0,0003 nD	0,0001 nD

Scope of application: Wine

The following models are particularly suitable for the measurement of the sugar content in fruit.

It indicates the expected °Alcohol of the fruit. The degree of ripeness of fruit (fruit-sugar) can also be determined, such as e.g. grapes. Alternatively the display can be switched to show Brix.

The main scope of applications is:

- Agriculture: Wine-growing (viticulture) and fruit-growing
- Wine-production
- Must and alcohol production



°Oe = Degree Oechsle, °KMW = Klosterneuburger Most Waage

Model	Scales	Measuring range	Accuracy	Division
KERN				
ORM 1WN	Oechsle	0 - 150 °Oe	± 2 °Oe	1 °Oe
	Percentage by volume	0 - 22 %	± 0,2 %	0,1 %
	KMW (Babo)	0 - 25 °KMW	± 0,2 °KMW	0,1 °KMW
	Brix	0 - 50 %	± 0,2 %	0,1 %
ORM 2WN	Oechsle France	0 - 230 °Oe	± 2 °Oe	1 °Oe
	Percentage by volume	0 - 22 %	± 0,2 %	0,1 %
	KMW (Babo)	0 - 25 °KMW	± 0,2 °KMW	0,1 °KMW
	Brix	0 - 50 %	± 0,2 %	0,1 %



Digital Refractometer KERN ORM-CO · ORM-UN

Scope of application: Coffee

The following models are particularly suitable for measuring the dissolved solids (TDS) in coffee to determine or compare the strength of a cup of coffee. For roasting plants, the TDS% value is used to determine the solubility level of a roast and to control the quality. Alternatively the display can be switched to show Brix or the refractive index.

The main scope of applications is:

- Coffee industry
- Coffee roasting plants
- Coffee competitions



Modell	Scales	Measuring range	Accuracy	Division
KERN				
ORM 1CO	Coffee TDS 1	0 – 25 %	± 0,2 %	0,1 %
	Brix	0 – 50 %	± 0,2 %	0,1 %
	Refractive index	1,3330 – 1,4200 nD	± 0,0003 nD	0,0001 nD
ORM 2CO	Coffee TDS 2	0,00 – 25,00 %	± 0,2 %	0,01 %
	Brix	0,00 – 30,00 %	± 0,2 %	0,01 %
	Refractive index	1,3330 – 1,4200 nD	± 0,0003 nD	0,0001 nD

Scope of application: Urine

The following models are particularly suitable for the measurement of the specific gravity (sg) in urine, the quantity of serum (serumproteine) in urine (doping control among athletes), and the refractive index.

The main scope of applications is:

- Hospitals
- Doctor's surgeries/Physicians
- Medical training institutions
- Nursing homes
- Sports medicine (doping test)
- Veterinary



Model	Scales	Measuring range	Accuracy	Division
KERN				
ORM 1UN	Urine (spec. gravity)	1,000 – 1,050	± 0,002	0,001
	Serum protein	0 – 12 g/100 ml	± 0,2 g/100 ml	0,1 g/100 ml
	Brix	0 – 50 %	± 0,2 %	0,1 %
	Refractive index	1,3330 – 1,4200 nD	± 0,0003 nD	0,0001 nD
ORM 2UN	Urine (s. g. dog)	1,000 – 1,060	± 0,002	0,001
	Urine (s. g. cat)	1,000 – 1,060	± 0,002	0,001
	Brix	0 – 50 %	± 0,2 %	0,1 %
	Refractive index	1,3330 – 1,4200 nD	± 0,0003 nD	0,0001 nD

Digital Refractometer KERN ORM-CA

Scope of application: Industry/Automotive

The following models are particularly suitable for the measurement and determination of AdBlue®, glycol concentration ethylene (EG) and propylene (PG), battery fluid (BF), urea, the freezing point of windscreens wash water (CW). Furthermore these models are suitable for the measurement of thermal exchange systems. Alternatively the display can be switched to show Brix or the refractive index.



Model	Scales	Measuring range	Accuracy	Division
KERN				
ORM 1CA	Wash water	(-60) - 0 °C	± 0,5 °C	0,1 °C
	AdBlue®	0 - 51 %	± 0,2 %	0,1 %
	Battery fluid	1,000 - 1,500 kg/l	± 0,005 kg/l	0,001 kg/l
	Brix	0 - 50 %	± 0,2 %	0,1 %
ORM 2CA	Refractive index	1,3330 - 1,4200 nD	± 0,0003 nD	0,0001 nD
	Ethylene glycol (%)	0 - 100 %	± 0,5 %	0,1 %
	Ethylene glycol (° C)	(-50) - 0 °C	± 0,5 °C	0,1 °C
	Propylene glycol (%)	0 - 100 %	± 0,5 %	0,1 %
	Propylene glycol (° C)	(-60) - 0 °C	± 0,5 °C	0,1 °C
	Brix	0 - 90 %	± 0,2 %	0,1 %



360° rotatable microscope head 360°	Fluorescence illumination for compound microscopes With 100W mercury lamp and filter	Integrated scale In the eyepiece	Battery operation Ready for battery operation. The battery type is specified for each device.
Monocular Microscope For the inspection with one eye	Fluorescence illumination for compound microscopes With 3 W LED illumination and filter	SD card For data storage	Battery operation rechargeable Prepared for a rechargeable battery operation
Binocular Microscope For the inspection with both eyes	Phase contrast unit For a higher contrast	USB 2.0 interface For data transmission	Plug-in power supply 230V/50Hz in standard version for EU. On request GB, AUS or USA version.
Trinocular Microscope For the inspection with both eyes and the additional option for the connection of a camera	Darkfield condenser/ unit For a higher contrast due to indirect illumination	USB 3.0 interface For data transmission	Integrated power supply unit Integrated in microscope. 230V/50Hz standard EU. More standards e.g. GB, AUS or USA on request.
Abbe Condenser With high numerical aperture for the concentration and the focusing of light	Polarising unit To polarise the light	WIFI data interface: For transmitting of the picture to a mobile display device	Package shipment The time required to manufacture the product internally is shown in days in the pictogram.
Halogen illumination For pictures bright and rich in contrast	Infinity system Infinity corrected optical system	HDMI digital camera For direct transmitting of the picture to a display device	Pallet shipment The time required to manufacture the product internally is shown in days in the pictogram.
LED illumination Cold, energy-saving and especially long-life illumination	Zoom magnification For stereomicroscopes	PC software To transfer the measurements from the device to a PC.	
Incident illumination For non-transparent objects	Auto-focus For automatic control of the focus level	Automatic temperature compensation For measurements between 10 °C and 30 °C	
Transmitting illumination For transparent objects	Parallel optical system For stereomicroscopes, enables fatigue-proof working	Protection against dust and water splashes IPxx: The type of protection is shown in the pictogram cf. DIN EN 60529:2000-09, IEC 60529:1989+A1:1999+A2:2013	
Fluorescence illumination For stereomicroscopes			

Abbreviations

C-Mount	Adapter for the connection of a camera to a trinocular microscope	SLR camera	Single-Lens Reflex camera
FPS	Frames per second	SWF	Super Wide Field (Field number at least Ø 23 mm for 10× eyepiece)
H(S)WF	High (Super) Wide Field (Eyepiece with high eye point for wearers of glasses)	W.D.	Working Distance
LWD	Long Working Distance	WF	Wide Field (Field number up to Ø 22 mm for 10× eyepiece)
N.A.	Numerical Aperture		